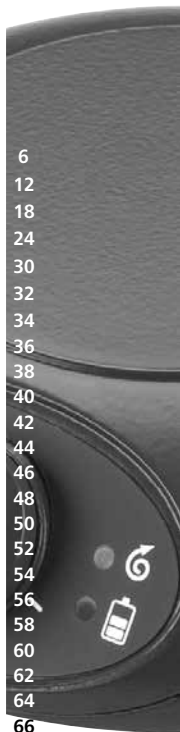
# IKA

designed for scientists

## IKA topolino mobil

Betriebsanleitung	DE	6
Operating instructions	EN	12
Mode d'emploi	FR	18
使用说明	ZH	24
Indicaciones de seguridad	ES	30
Veiligheidsinstructies	NL	32
Norme di sicurezza	IT	34
Säkerhetsanvisningar	SV	36
Sikkerhedshenvisninger	DA	38
Sikkerhedsanvisninger	NO	40
Turvallisuusohjeet	FI	42
Normas de segurança	PT	44
Wskazówki bezpieczeństwa	PL	46
Bepečnostní upozornění	CS	48
Biztonsági utasítások	HU	50
Varnostna navodila	SL	52
Bezpečnostné pokyny	SK	54
Ohutusjuhised	ET	56
Drošības norādes	LV	58
Saugos reikalavima	LT	60
Указания за безопасност	BG	62
Instrucțiuni de securitate	RO	64
Υποδείξεις ασφαλείας	EL	66

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## Contents

	Page
Safety instructions	12
Correct use	14
Unpacking	14
Commissioning	14
Operating	15
Accessories	16
Maintenance and cleaning	16
Warranty	17
Technical data	17

## Safety instructions

### • Read the operating instructions in full before starting up and follow the safety instructions.

Keep the operating instructions in a place where they can be accessed by everyone.

- Ensure that only trained staff work with the appliance.
- Use only the AC adapter supplied.
- Observe the following safety instructions for the handling of batteries. Considerable material and health damage and personal injury can be caused by:
  - Leaking or exploding batteries
  - Escaping of the toxic or corrosive substances contained in batteries
  - Use only the rechargeable battery types recommended in the technical data in the appliance!  
Do not use other battery types and on no account non-rechargeable batteries.
  - Never mix different models, types or capacities of

batteries.

- Be sure to observe the polarity indicated in the battery compartments! The side of the battery with the plus sign (+) must be aligned with the correspondingly marked side of the battery compartment.
- Do not charge batteries that have leaked or that are discoloured, deformed or damaged in any other way.
- Leaking or damaged batteries must be disposed of while observing the normal precautionary measure (protective gloves).
- Do not open, dismantle, pierce, modify or throw the batteries and do not expose them to unnecessary blows.
- Do not expose the batteries to fire or temperatures above 50 °C. Do not store the batteries in hot places (e.g. in the car in direct sunlight).
- Do not charge the batteries when the ambient temperature is below 0 °C or above 40 °C.
- Do not immerse the batteries in water or other liquids.
- Objects made of metal or other conductive materials (e.g. keys, coins, tools, rings) can cause a short-circuit in the batteries. Never store the batteries together with such objects.
- Always replace the batteries as a complete set. Never use old and new batteries together.
- Never connect the batteries directly to a power source (e.g. a domestic plug socket or a car battery).
- Never place the batteries in a microwave oven or in pressurised or vacuum vessels.
- If you charge the batteries in an external charger, ensure that the admissible charging current for these

- batteries is not exceeded
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- **Caution - Magnetism!** Effects of the magnetic field have to be taken into account (e.g. data storage media, cardiac pacemakers...).
- Wear your personal protective equipment in accordance with the hazard category of the media to be processed. Otherwise there is a risk from:
  - splashing and evaporation of liquids
  - Escape of toxic or inflammable gases
- The appliance is not suitable for manual operation.
- Set up the appliance in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- The feet of the appliance must be clean and undamaged.
- Position the knob at the left stop before starting up. Gradually increase the speed. Reduce the speed if
  - the medium splashes out of the vessel because the speed is too high
  - the appliance is not running smoothly
  - the container moves on the base plate.
- Check the appliance and accessories beforehand for damage.
- Do not use damaged components. Only replace damaged parts with spare parts identical to the original in function and quality.
- Beware of and avoid hazards due to
  - flammable materials
  - glass breakage
  - incorrect container size
  - overfilling of media
  - unsafe condition of container or of the complete assembly.

- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any energy produced in other ways, e.g. through light irradiation.
- It may be possible for wear debris from rotating accessory parts to reach the material being processed.
- Do not use the device in explosive atmospheres, it is not EX-protected.
- With substances capable of forming an explosive mixture, appropriate safety measures must be applied, e.g. working under a fume hood.
- To avoid body injury and property damage, observe the relevant safety and accident prevention measures when processing hazardous materials.
- The appliance must only be disconnected from the mains supply by pulling out the plug-in power supply unit.
- Safe operation is guaranteed only with the use of original IKA accessories.
- The device will automatically restart following any interruption to the power supply.
- The appliance may heat up when in use.
- The voltage stated on the type plate must correspond to the mains voltage.
- Protect the appliance and accessories from bumps and impacts.
- The appliance and the power supply unit are maintenance-free and must not be opened.

- Dispose of old batteries and defective appliances in accordance with the statutory regulations.

## Correct use

### • Use

- For mixing of liquids independently of a mains power supply

### • Range of use

Indoor environments similar to that a laboratory of research, teaching, trade or industry area.

## Unpacking

### • Unpacking

- Please unpack the device carefully
- In the case of any damage a detailed report must be set immediately (post, rail or forwarder).

### • Contents of package

- Magnetic stirrer
- Plug-in power supply unit
- Four batteries (AA)
- Operating instructions

## Commissioning

- Set the knob to the left stop (you will hear and feel it clicking into place)

### Insert the batteries

- Disconnect the AC adapter from the appliance before inserting the batteries.
- The battery compartments are located on the underside of the appliance. Remove the battery compartment cover by pressing the two tabs on the side.
- Pay attention to the correct polarity when inserting the batteries. The polarity is indicated in the battery compartment.

- Fit the battery compartment cover again and press to close. The cover engages audibly.

### Caution!

Use only the rechargeable Type AA batteries recommended by the manufacturer. Never use conventional, non-rechargeable batteries as considerable material and health damage and personal injury cannot be ruled out. The batteries could explode or be destroyed. Furthermore, the appliance may be damaged or destroyed. Use only four rechargeable batteries of the same type and same capacity.

### Connect the AC adapter

- Insert the round plug of the AC adapter to the socket on the rear of the appliance
- Connect the AC adapter to a plug socket

The batteries are charged as soon as the AC adapter is plugged in.

### Charging the batteries

You can continue to use the appliance stirrer drive during the entire charging process.

The charging cycle is performed automatically in the following steps:

- Step 1: Rapid charging - red charge indicator  flashing  
The duration of the charging cycle depends on the discharge state and capacity of the battery and can be up to 3 hours. Charging of the batteries during the cycle: approx. 0 - 80%.
- Step 2: Normal charging - red charge indicator  steady light

The duration of this charging cycle depends on the discharge state and capacity of the battery and can be up to 30 minutes. In this state the appliance can already be used independently of a mains power supply.

Charging of the batteries during the cycle:

approx. 80 - 100%.

- Step 3: Trickle charge - red charge indicator  not lit  
The batteries are completely charged, the charge is maintained. The appliance can be used independently of a mains power supply. Charging of the batteries during the cycle: approx. 100%.

This cycle starts after every disconnection from the mains power supply!

**During initial use of the appliance or the first use of a new set of batteries:** New, uncharged batteries have only a precharge and must be charged for at least 4 hours before using the appliance for the first time. In order to achieve the maximum capacity, the first charge cycle of the batteries should not be interrupted!

Batteries become warm during charging; this is a normal phenomenon and is not dangerous. Batteries are subject to self-discharging. If the appliance is seldom used, the batteries should nevertheless be charged roughly 1x per month to prevent a deep discharge.

Do not leave the batteries or the appliance unused or uncharged for prolonged periods. If batteries are not used for longer than a few months, they may become unserviceable and have to be replaced. If the appliance is not to be used for a prolonged period, remove the batteries and store them in a dry place at normal room temperature.

Deep discharging of the batteries should always be avoided as this negatively affects the service life of the batteries. If the speed of the appliance drops noticeably, the batteries should be recharged.

Batteries are wear parts and are subject to ageing during normal use. A reduction in the capacity and hence in the useful operating time during the course of time is normal and not a fault. The warranty granted on the appliance therefore does not apply to the batteries. In order to maintain the full running time of the appliance, batteries should be replaced after approx. 500 - 600 charge cycles or 2 years.

The appliance can be operated with rapid-charging NiCad or NiMH AA batteries with capacities from 1000 mAh to 2700 mAh, but always only with four identical batteries at the same time. NiMH batteries are to be recommended due to their higher capacity. The charging and operating times shown are indicative values and related to the batteries supplied with the appliance.

## Operation

The appliance can be operated permanently with the AC adapter or with batteries independently of a mains power supply. If the appliance is operated with an AC adapter, the batteries automatically take over the power supply in the event of a mains power failure. During battery operation, the operating time independently of a mains power supply depends on the condition of the batteries. If all the batteries are intact and fully charged, operating times of > 8 hours are achieved with the batteries supplied with the appliance.

- Check before use whether the charge state of the batteries is sufficient or whether the AC adapter is plugged in.
- Pour the liquid to be stirred into a suitable vessel and insert an appropriate magnetic bar. The vessel must be permeable to a magnetic field.
- Position the vessel in the middle of the set-up plate.
- Switch on the motor at the rotary knob, the green operating indicator  lights up continuously.
- Set the speed with the rotary knob. Turning to the right increases the speed, turning to the left decreases the speed.
- After stirring, turn the rotary knob completely to the left until it engages audibly and perceptibly. In the left-hand end position, the motor is switched off and the green operating indicator is out.
- If the appliance does not function or exhibits inadequate performance, please check whether
  - All the batteries are inserted with the correct polarity,
  - The batteries are sufficiently charged.

## Accessories

- Magnetic stirring bars: max.  $\varnothing$ 6mm x 30mm
- Recommended batteries:
  - Mignon rapid-charging (AA) batteries of types NiCad or NiMh with 1.2 V rated voltage and a capacity from 1000 mAh to 2700 mAh.

Spare parts list see [www.ika.com](http://www.ika.com)

## Maintenance and cleaning

The device is maintenance-free

Only the batteries are wear parts and must be replaced when necessary. The batteries are worn out when despite a complete charge, only short operating times are achieved. Always replace all 4 batteries at the same time.

### Cleaning

- For cleaning disconnect the mains plug!
- Use only cleaning agents which have been approved by IKA to clean IKA devices.
  - These are water (with tenside) and isopropanol.
- Wear protective gloves during cleaning the devices.
- Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.
- Do not allow moisture to get into the device when cleaning.
- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.

## Warranty

In accordance with IKA warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine directly to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

## Technical data

Rated voltage	<b>VAC</b>	100 - 240
Rated frequency	<b>Hz</b>	50/60
Absorbed power	<b>W</b>	10
Power input	<b>W</b>	7.2
Batteries		4x Type Mignon (AA), NiCad or NiMh, 1.2 V, 1 Ah - 2.7 Ah
Speed (adjustable)	<b>rpm</b>	0 / 500...1800 $\pm$ 25%
Operating time	<b>%</b>	100
Ambient temperature	<b>°C</b>	+5 to +40
Ambient humidity (relative)	<b>%</b>	80
EN 60529 protection class		IP 21
Pollution degree		2
Overvoltage category		II
Protection class		II
Transport temperature	<b>°C</b>	-25 to +50
Storage temperature	<b>°C</b>	-10 to +50
Operation at a terrestrial altitude	<b>m</b>	max. 2000 above sea level
Set-up plate	<b>mm</b>	$\varnothing$ 80
Stirring bar (max.)	<b>mm</b>	$\varnothing$ 6 x 30
Max. stirring quantity	<b>ml</b>	250
Dimensions (D x H)	<b>mm</b>	140 x 42
Weight	<b>kg</b>	0,6 (incl. battery and AC adapter)

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**IKA Vietnam Company Limited**
**IKA India Private Limited**